

# Kimtech Pure\* A5 Sterile Cleanroom Apparel Gowning Procedure

## Before Gowning

### Step 1

(Pre-Entry) Don Hair Net and Shoe Covers after removing all jewelry and cosmetics.

### Step 2

(Gowning) Wash hands and gown first pair of sterile gloves. Sanitize gloves after gowning each article if required.

### Step 3

Apply mask and hood assuring a snug fit.

### Step 4

Open vacuumed-packed apparel. Tear at notched edge.

### Step 5

Grasp the blue line. Located on the inside middle back.



### Step 6

Gently unfold coverall utilizing blue indicator line. Arms and legs are pre-drawn and snapped in place. Garment is already folded inside-out and unzipped.



*Snaps allow gathered-up arms and legs to expand during gowning.*

## Begin Gowning

### Step 7

Hold garment at waist.



### Step 8

Put one leg in and point toe through opening until snap releases.



### Step 9

Do the same with the other leg.



### Step 10

Insert one arm and extend until snap releases.



### Step 11

Do the same with the other arm.



### Step 12

Slip thumbs through thumb loops.



### Final Step

Cross legs and zip up coverall.



### After A5 Sterile Cleanroom Apparel is gowned

- Add boot covers.
- Complete gowning by adding goggles and a second pair of sterile gloves.



Exceptional Workplaces<sup>®</sup>

For more information, contact your local Kimberly-Clark Professional<sup>®</sup> Sales Representative at 1-800-255-6401 or visit [www.kimtech.com](http://www.kimtech.com).

